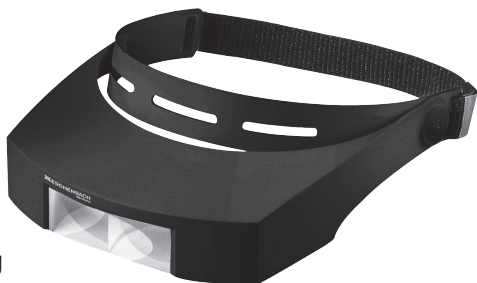
laboCOMFORT

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User Manual

Mode d'emploi

Istruzioni d'uso

Manual de instrucciones

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Bruksanvisning

Käyttöopas

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English

Congratulations on purchasing this ESCHENBACH headband magnifier - a quality product 'Made in Germany'.

Intended use

laboCOMFORT is used where precise work of detection of the smallest details is required. Depending on the field of application, lens parts of various magnifications can be fitted and different working distances (distance lens to object) can be applied.

Safety information

▶ **Danger of fire!**

Lenses in optical devices can cause significant damage by focusing light to generate heat if handled or stored improperly. Be careful never to leave optical lenses lying in direct sunlight!

▶ **Danger of glare and injury!**

Never look directly into the sun or other bright light sources using optical devices!

▶ **Also inform other people and children especially of this safety measure!**

- ▶ **laboCOMFORT may not be used while driving automobiles.**
- ▶ **Protect your laboCOMFORT from impacts and excessive heat. Never lay your laboCOMFORT on a radiator or in direct sunlight.**
- ▶ **The product is a Class I medical device (non-sterile, without measuring function) and complies with Regulation (EU) 2017/745. For use of the product outside the EU member states, the respective country-specific directives must be observed.**

Operating elements

- 1 Lens part
- 2 Magnifier case
- 3 Forehead support with air vents
- 4 Elastic strap with Velcro fastener

Adjusting the headband magnifier

The strap can be easily adapted to any head size by adjusting the Velcro fasteners on each side of the strap. The forehead support has gaps so that air can circulate. The magnifier case of the headband magnifier can be set to the most comfortable position through various snap-in positions. This allows for a comfortable head position. When the headband magnifier is not in use it does not need to be taken off completely but can be simply swiveled up.

Exchanging the lens

To exchange the lens pull it down out of the frame of the magnifier case. A different lens can then be snapped into the frame. Just make sure that the curved surface of the lens is on the side away from the eye. The magnification of the lens is stamped into the lens' adapter. The headband magnifier is available with various magnifications. There is a certain working distance (distance between lens and viewed object) for each magnification that should be kept to get an optimal image quality. Extra individual lenses can be ordered by the following order numbers:

Order No.	Magnification	Working Distance
16455	1.7 ×	Approx. 400 mm
16451	2.0 ×	Approx. 250 mm
16452	2.5 ×	Approx. 180 mm
16453	3.0 ×	Approx. 160 mm

Care instructions

All ESCHENBACH lenses are manufactured from PXM[®], an optically pure plastic with user-friendly characteristics:

- unbreakable
- lighter than glass
- optically pure for best viewing quality

PXM[®] lenses are coated in a special process, are scratch-resistant, slightly bloomed, and given a permanent anti-static coating.

Apparent hairline cracks that may be visible along the lens' edge do not impair the optical image. They are flow lines of the material that occur during production.

- ▶ Do not use alcohol or organic solvents to clean headband magnifiers.
- ▶ Do not use any soap solutions containing softeners or any abrasive cleaning products. These may ruin the lenses.
- ▶ Clean the lenses with a soft cotton or linen cloth (such as cleaning cloths for eyeglasses). Moisten the cleaning cloth with lukewarm water to remove more persistent stains.

Warranty

Within the framework of the applicable statutory provisions, we provide a warranty for the functioning of the product described in this manual with

regard to faults arising in connection with manufacturing errors or material errors. In event of damage due to improper handling, including damage from a fall or impact, no warranty claim exists. Warranty claims are only accepted upon presentation of the sale receipt!

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Eschenbach Optik GmbH